

# Ginsenoside Analysis in Ginseng extract

## 1. Analytical Condition

### YL9100 Series HPLC

Mobile Phase : ACN:Water=Gradient

Flow Rate : 1.0ml/min

Column Oven :35°C

Detector : UV/Vis 203nm

Column : C18 (4.6\*150mm, 3.5um)

Injection Volumn : 20ul sample loop

Gradient Program	Water	ACN
Initial	80	20
10.00	80	20
40.00	68	32
48.00	58	42
50.00	0	100
60.00	0	100
62.00	80	20

## 2. Chromatogram

